# **INSTRUCTION MANUAL**





# HI6321 Advanced Conductivity Benchtop Meter

Hanna Instruments Inc., 584 Park East Drive, Woonsocket, RI 02895 USA www.hannainst.com

# Dear Customer,

Thank you for choosing a Hanna Instruments® product.

Please read this instruction manual carefully before using this instrument as it provides the necessary information for correct use of this instrument, as well as a precise idea of its versatility.

If you need additional technical information, do not hesitate to e-mail us at tech@hannainst.com. Visit www.hannainst.com for more information about Hanna Instruments and our products.

# **TABLE OF CONTENTS**

1.	Prelim	inary Examination	. 3
2.	Safety	Measures	. 3
3.	User li	nterface — Icons	. 4
4.	Genero	al Description & Intended Use	. 6
5.	Specifi	cations	. 8
	5.1.	Instrument	. 8
	5.2.	HI7631233 Probe	11
6.	Function	onal & LCD Description	12
	6.1.	Main Unit	12
	6.2.	Probe	14
7.	Getting	g Started	15
	7.1.	Installing the Holder	15
	7.2.	Using the Holder	16
	7.3.	Connecting the Probe, Keyboard, Printer	16
	7.4.	Powering the Unit & Selecting Operating	
		Language and Regional Preferences	
	7.5.	Basic Operations	17
8.	Systen	n Menu Items	18
		Users	
	8.2.	System Settings	22
		Measurement Settings	
	8.4.	Log Recall	30
	8.5.	Help	36
9.	Measu	rement & Probe Setup Menu	37
	9.1.	Calibration	37
	9.2.	Reading	38

9.3. Temperature	40
9.4. View	42
9.5. Alarms	45
9.6. Logging	46
9.7. Profiles	47
10. Calibration	48
10.1. Conductivity Calibration	48
10.2. Procedure	
10.3. Salinity Calibration	50
11. Measurement	51
11.1. Measurement Tips	52
11.2. Direct Readings	52
11.3. Direct / Autohold Readings	52
12. Logging	53
12.1. Automatic Logging	
12.2. Manual Logging	54
12.3. Autohold Logging	55
13. Maintenance	56
14. Error Messages	57
15. Software Update	58
16. Accessories	
Certification	59
Recommendations for Users	59
Warranty	59
Regulatory Notices for the Wi-Fi Module	60

All rights are reserved. Reproduction in whole or in part is prohibited without the written consent of the copyright owner, Hanna Instruments Inc., Woonsocket, Rhode Island, 02895, USA. Hanna Instruments reserves the right to modify the design, construction, or appearance of its products without advance notice.

# **1. PRELIMINARY EXAMINATION**

HI6321 is Hanna's<sup>®</sup> advanced benchtop conductivity meter with a large touch screen display and streamlined design. Each HI6321 is supplied with:

- HI7631233 EC and resistivity probe
- HI764060 electrode holder with following accessories:
  - base plate (with integrated pivot pin) and screw (requires installation)
  - cable holder clip, attached
  - electrode holder with adapter, attached
- EC calibration starter kit consisting of:
  - $\circ$  1413  $\mu$ S/cm standard solution (4 sachets)
  - $\circ$  12880  $\mu$ S/cm standard solution (2 sachets)
  - $\circ$  5000  $\mu$ S/cm standard solution (2 sachets)
- Electrode rinse solution (2 sachets)
- 24 VDC power adapter
- USB-C to USB-A cable
- Probe quality certificate
- Quick reference guide
- Instrument quality certificate

**Note:** Save all packing material until you are sure the instrument works correctly. Any damaged or defective item must be returned in its original packing material with the supplied accessories.

# 2. SAFETY MEASURES

# Handling & Usage Precautions

The unit, while not fragile, can be damaged by improper handling and usage.

- Transport the unit with all cables removed.
- Keep the unit on a stable and even surface, away from contact with liquid.
- Avoid excessive dirt and dust.
- Protect the unit from contact with food, oils, and chemicals.
- If the device becomes wet, gently wipe the exterior with a clean, dry cloth.
- Keep away from direct sunlight.
- Use in a safe place that is appropriate to application requirements.
- Use attachments and accessories specified in this manual only.
- Capacitive touchscreen and buttons operate without applying pressure.
- Do not puncture the capacitive touchscreen or drop the unit.
- Do not use the device near heat sources.
- Do not place objects on top of the device.
- Do not insert objects into the ports, spaces around keys, other than the intended cable, USB drive.

# 3. USER INTERFACE - ICONS

Capacitive keys	Description
<	Back — return to a previous hierarchical menu level
0	Home – access to measurement screen & configured profile
=	<b>Menu</b> — access to Users, System Settings, Measurement Settings, Log Recall, Help
Main menu	Description
<u>.</u>	Users — login & rights configuration
<b>e</b>	eq:System System System System Configuration, connectivity & printing items
Ô	Measurement Settings
$\bigcirc$	Log Recall — access logged measurement data
2	Help — access support
Measurement	Description
<b>\$</b>	Measurement Menu, accessed from measurement screen
	EC and resistivity probe
A	Autohold applied
EC / Res / TDS / Sal	Autohold, pending stable measurement
A / A	Warning on standby / active function
Logging	Description
# 00002 00:00:12	Start/Stop logging (current index and time since log start)
<b>1 1 1 1 1 1 1 1 1 1</b>	Manual logging (current index)
0	Triggers log session, pending next stable measurement
A	Autohold logging in progress
) ( <b>D</b>	Annotated text/Annotated text in use
0% 25% 50%	Available file storage capacity (at full capacity the icon is displayed blinking)

Log recall		Description		
	🔲 / 🗐	Table view — function active/not selected		
	🗠 / 🗠	Graph view — function active/not selected		
	<b>()</b> / <b>()</b>	Information view — function active/not selected		
Genero	ıl	Description		
	ð	Measurement profile		
	X	Background operation in progress		
	Ø	Active alarm set		
Un	stable Stable Autohold	Stability/Autohold indicator		
Connec	tivity & Printing	Description		
+	品	Connection established (tap for IP address)		
Ethernet	윰 0.5 sec 윰	Connection in progress		
· تن	Z	Connection error		
	<b>(</b> [r	Connection established (tap for IP address)		
Wi-Fi	🛜 0.5 sec 🛜	Connection in progress		
	All Contractions	Connection error		
ISB	B	USB-A or USB-C flash drive plugged in		
S		High-power consumption with the flash drive plugged in		
PC		PC connection established through USB-C port		
	¢	Printer connected — printing manual log option <b>enabled</b>		
Printer	Ē	Printer connected — printing manual log option <b>disabled</b>		
	<u><u></u></u>	Printer not recognized or no longer connected		

# 4. GENERAL DESCRIPTION & INTENDED USE

HI6321 is an advanced benchtop meter with a capacitive display, comprised of a housing and an integrated conductivity measurement module. Compact and easy to operate, the meter is delivered with HI7631233 EC and resistivity probe.

HI7631233 is a platinum, four-ring probe with an integral temperature sensor that adjusts the measured conductivity to a reference temperature by applying compensation algorithms:

- linear: when the temperature coefficient of variation has the same value for all measurement temperatures
- standard: for high-purity water measurements (  $>\!1\,M\Omega$  cm resistivity) and documented in ASTM Standard D5391-14

• **natural**: for natural ground, well, or surface water (water with similar composition) as per ISO7888 standard The result is reliable electrolytic conductivity, resistivity, TDS (Total Dissolved Solids), or Salinity (expressed in percent, PSU, or ppt units).

**TDS** is a calculated value based on the conductivity of the solution (TDS = factor  $\times$  EC<sub>25</sub>). A TDS factor is a conversion factor used to change an EC measurement to a ppm measurement.

**Sal PSU**: The practical salinity of seawater relates the ratio of electrical conductivity of a normal seawater sample at 15 °C and 1 atmosphere to a potassium chloride solution (KCI) with a mass of 32.4356 g/Kg water at the same temperature and pressure. Under these conditions the ratio is equal to 1 and S = 35. The practical salinity scale may be applied to values 0 through 42.00 PSU at temperatures between 0 to 35 °C.

**Sal ppt**: measurements expressed in ppt are based on the Natural Seawater Scale that extends from 0.00 to 80.00 g/L and covers 10 to 31 °C temperature range. It determines the salinity based upon a conductivity ratio of sample to standard seawater at 15 °C and an approximate salinity value of 35 in seawater.

Sal %: in this scale 100 % salinity is equivalent to roughly 10 % solids.

This system responds to a complex range of measurement and monitoring requirements, providing accuracy, reproducibility, and reliability.

HI6321 is supplied with an electrode holder that has a flexible arm. The holder can be mounted quickly and provides secure support for electrodes while taking measurements in sample containers.

This benchtop meter supports:

- USB type A support for USB stick, printer, keyboard
- USB type C support for USB stick, PC connection

The user can select between five different views.

- Basic measurement configuration
- Simple GLP with calibration information
- Full GLP with probe status and calibration point details
- Live updated, interactive graph
- Tabulated data with date, time, and notes

# Capacitive touch screen with multi-touch support

The benchtop unit has a 7-inch color display with resolution of  $800 \times 480$ . The capacitive, multi-touch screen supports video playback and data plotting.

# **Main Features**

## **Calibration & Measurement**

- Measure
  - Conductivity  $-\mu$  S/cm or mS/cm
  - Resistivity Ω·cm, kΩ·cm, MΩ·cm
  - ► TDS ppm or ppt
  - ▶ Salinity ppt, PSU, %
- Application-specific profiles for quick and direct measurement without the need to update the sensor and system settings
- Active log during measurement
- Measurement stability indicator (using the Stability Criteria setting)
- Reading modes: direct and direct/autohold
- Temperature source (Automatic or set manually)
- Audible and/or alarm messages for measurements outside threshold limits
- Galvanic isolation for conductivity measurement module
- Standard (up to)**four-point conductivity** calibration with standards for cell factor determination and 0  $\mu$ S/ cm for offset:
  - 84  $\mu$ S/cm, 1413  $\mu$ S/cm, 5.0 mS/cm,12.88 mS/cm for the 0.1/cm cell
  - 80.0 mS/cm, 111.8 mS/cm additional standards for the 1.0/cm cell
- Standard single point salinity calibration in 100 % salinity standard, with Salinity scale set to %
- Non-volatile memory saves data and settings

# Logging

- Data log collection of at least 1 000 000 data points (with time and date stamp)
- Logging types: manual, automatic, autohold
- Sample ID for manual and Autohold data

# Connectivity features & services

- Transfer logged data to a USB thumb drive
- Log files that include measurements and calibration data (as .CSV file)
- FTP and email for log export via Ethernet and Wi-Fi connection
- Download logs using the benchtop's embedded web server
- USB type A for USB stick, printer, keyboard
- USB type C for USB stick and PC connection

# User-support feature

• Help section – brief overview of instrument's main functionalities and features

# 5. SPECIFICATIONS

# 5.1. INSTRUMENT

		0.000 to 9.999 µS/cm
		10.00 to 99.99 µS/cm
	Range	100.0 to 999.9 µS/cm
	Kuliyo	1.000 to 9.999 mS/cm
		10.00 to 99.99 mS/cm
		100.0 to 1000.0 mS/cm
		0.001 µS/cm
		0.01 µS/cm
	Resolution	0.1 µS/cm
	Kesolulion	0.001 mS/cm
		0.01 mS/cm
		0.1 mS/cm
	Accuracy	$\pm 1$ % of reading or $\pm 0.010 \mu$ S/cm, whichever is greater
	Cell constant	0.0500 to 200.0000 /cm
		Automatic
	Calibration type	Manual
	Calibration points	Single
Conductivity		Multiple
	Calibration solutions	84.00 µS/cm
		1.413 mS/cm
		5.000 mS/cm
		12.88 mS/cm
		80.00 mS/cm
		111.8 mS/cm
		Disabled
	Calibration reminder	Daily: 0 min. to 23 hours and 59 min.
		Periodic: 1 min. to 30 days, 23 hours and 59 min.
		Linear
	Temperature	Natural
	compensation	Standard
	·	Disabled
	Reference	Range 5.0 to 30.0 °C (41.0 to 86.0 °F)
	temperature	Resolution 0.1 °C/0.1 °F
	Temperature coefficient	0.00 to 10.00 %/°C

		1.0 to 99.9 Ω•cm				
		100 to 999 Q • cm				
		1.00 to 9.99 KΩ•cm				
	Range	10.0 to 99.9 KΩ • cm				
		100 to 999 KΩ•cm				
		1.00 to 9.99 MQ • cm				
		10.0 to 100.0 MΩ • cm				
Resistivity		0.1 Ω•cm				
		1 Ω•cm				
		0.01 KΩ•cm				
	Resolution	0.1 KΩ•cm				
		1 KΩ•cm				
		0.01 MΩ•cm				
		0.1 MΩ•cm				
	Accuracy	$\pm 1$ % of reading or $\pm 1$ $\Omega$ · cm, whichever is greater				
		0.000 to 9.999 ppm				
		10.00 to 99.99 ppm				
		100.0 to 999.9 ppm				
	Range	1.000 to 9.999 ppt				
		10.00 to 99.99 ppt				
Total		100.0 to 400.0 ppt				
Dissolved		actual TDS (with 1.00 factor)				
Solids		0.001 ppm				
(TDS)		0.01 ppm				
	Resolution	0.1 ppm				
	Resolution	0.001 ppt				
		0.01 ppt				
		0.1 ppt				
	Accuracy	$\pm 1$ % of reading or $\pm 0.01$ ppm, whichever is greater				
		0.00 to 42.00 PSU - Practical Scale				
	Range	0.00 to 80.00 ppt - Natural Sea Water				
	-	0.0 to 400.0 % - Percent Scale				
Salinity	Develution	0.01 for Practical Scale / Natural Sea Water				
	Resolution	0.1 % for Percent Scale				
	Accuracy	$\pm 1$ % of reading				
	Calibration	1 point for percent scale, using 100 % salinity calibration solution				

	<u> </u>				
	Range	-20.0 to 120.0 °C/ -4.0 to 248.0 °F/ 253.2 to 393.2 K			
Temperature	Resolution	0.1 °C/0.1 °F/0.1 K			
	Accuracy	$\pm 0.2 \text{ °C} / \pm 0.4 \text{ °F} / \pm 0.2 \text{ K}$			
	Calibration	Single point, adjustable			
	Modes	Direct Direct/Autohold			
Reading	Stability criteria	Accurate Medium Fast			
	Basic	Conductivity, Resistivity, TDS, Salinity, Temperature measurement data Measurement profile (if enabled) Stability status			
Views	Simple GLP	Basic view information Last calibration date and offset			
	Full GLP Conductivity & Salinity	Simple GLP information and calibration point details Cell constant, reference and source temperature			
	Graph (Plot)	Measurement versus time graph (panned or zoomed)			
	Table	Measurements updated every second (displayed in table)			
	Туре	Automatic, Manual, Autohold			
	Number of records	50 000 maximum per file Stores at least 1 000 000 data points per user			
Logging	Automatic interval	1, 2, 5, 10, 30 seconds 1, 2, 5, 10, 15, 30, 60, 120, 150, 180 minutes			
	Sample ID	Incremental mode or manual			
	Export option	.CSV file format			
Users		Up to 9 users and the default administrator account			
	USB-A	2 ports for keyboard and/or printer input or USB thumb drive			
	USB-C	1 port for PC connectivity, USB-C type thumb drive and printer			
Connectivity	Wi-Fi & Ethernet	FTP Web server Log transfer and download Email			
	RS232	Connecting peripherals			
Power supply		DC adapter 100-240 VAC to 24 VDC 2 A			
Environment		0 -50 °C / 32 - 122 °F / 273 - 323 K Maximum 95% RH non-condensing			
Dimensions		205×160×77 mm (8.0×6.2×3.0")			
Weight		Approximately 1.2 kg (26.5 lbs.)			


# 5.2. HI7631233 PROBE

Range	0 to 1000 mS/cm		
Recommended operating temperature	—5 to 100 °C (23 to 212 °F)		
Temperature sensor	Built-in		
Cell constant	1 ± 15 %		
Cell type	Four-ring, platinum on glass		
Body	Polyetherimide (PEI)		
Wetted parts	Sensor body PVDF O-ring NBR		
Sensor diameter	Ø 12.0 mm		
Connection	DIN		
Cable length	1 m (3.3′)		


# 6. FUNCTIONAL & LCD DESCRIPTION

#### 6.1. MAIN UNIT

#### **Front View**

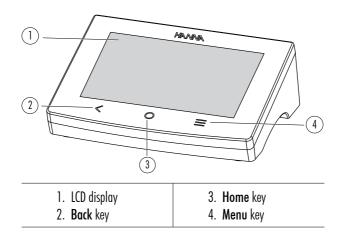

#### **LCD Description**

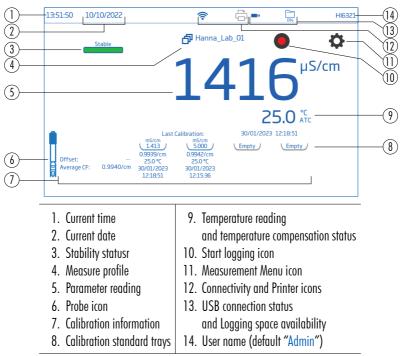

# Status Area

10:07:11 03/03/2022 🛜 📼 🛱 Admin


Status area runs horizontally across the top of the LCD screen and is continuously displayed after powering the unit.

- Left corner current time & date
- Middle connectivity status
- Right corner storage space availability & username

#### **Direct Keys**

lcon	Name	Function
<	Back	<ul> <li>returns the user to previous hierarchical menu level</li> <li>exit or escape function</li> </ul>
0	Home	<ul><li> access to measurement screen</li><li> exit or escape function</li></ul>
Ξ	Menu	• access to Users, System Settings, Measurement Settings, Log Recall, Help


# **Rear View**

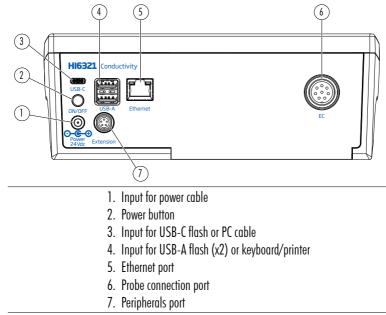

# 6.2. PROBE

The HI7631233 four-ring conductivity probe provides stable measurements over a wide measure range and does not require frequent calibrations.

An integral temperature sensor measures temperature and adjusts the measured conductivity to a reference temperature by applying compensation standards.

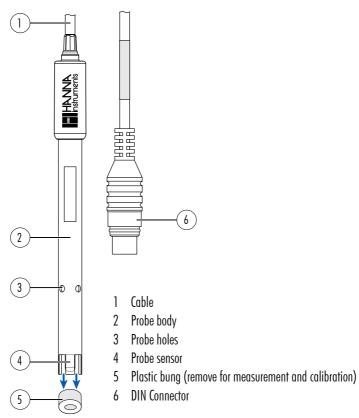


# 7. GETTING STARTED

# 7.1. INSTALLING THE HOLDER

# Attaching the base plate


- Take the HI764060 holder from the box.
- Identify the metal base plate (4) with the integrated pivot pin (5) and the screw (2).
- The plate may be attached to either side of the meter, left (L) or right (R).
- Place the meter face down on a clean, dry surface.
- Align the hole on the base plate (3) over the rubber foot (6R or 6L). The pivot pin (5) should be facing downward.
- Use a screwdriver (1) to tighten the screw (2) and attach the base plate to the meter.
- Position the meter with the display facing up.
- Slide the holder (7) over the pivot pin (5). A "slide in" motion is required to lock the arm into position.
- For increased arm rigidity, tighten the metal knobs on both sides of the holder.

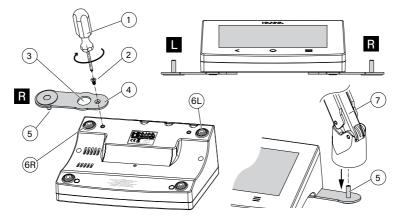

# Cable holder clip

The holder is delivered with a cable clip (attached) that secures several cables whilst allowing them to move freely with the arm motion.

- 1. To open the latch, press the clip inward while pulling up the latch.
- 2. To close the latch, lower latch over cable and snap closed. The latch snaps in position and secures the cables inside.

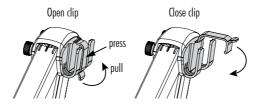

# **Getting Started**

To reattach the clip onto the electrode arm:

- 1. Align the clip's dovetail over the slot.
- 2. Gently push down to slide in position.

## Using the adapter

The arm holder is ended with a probe holder fitted with an adapter that has three different-sized apertures:

- center-front (reference temperature probe only)
- center-back (conductivity probe, with adapter only)
- left and right (conductivity probe, no adapter)

# 7.2. USING THE HOLDER

Use the holder for probe support and easy movement in and out of beakers and containers during calibration and sample measurement.

# 7.3. CONNECTING THE PROBE, KEYBOARD, PRINTER

# Connecting the probe

The HI7631233 probe is connected to the benchtop through a DIN connector, which makes attaching and removing the probe an easy process. With the meter off:

- Connect the probe to the DIN connection port, on the back of the meter. Align the pins and key then push the plug into the socket.
- Place the probe into the holder and secure the cable in the cable holder clip.

Note: Connector and plug must be connected firmly for the system to work correctly.

# Connecting a USB-A keyboard

Connect a USB's keyboard plug into the USB-A input on the back of the unit. Once connected the keyboard is automatically detected.

Users can attach a keyboard to set up calibration password, sample ID, user ID, or company/lab name.

# Connecting a printer

Hanna<sup>®</sup> aims to ensure meter compatibility with USB printers but can not ensure compatibility with all models. HI6321 can print directly to certain models of USB-dedicated printers with PCL printer language capability.

#### Printer components and requirements

- Printer, PCL driver compatible
- Cable
  - ▶ power cable
  - USB connector cable with two ends: type B connector (plugs into printer) type A connector (plugs into the USB port on the meter)



Side positioning

Center-back positioning




# 7.4. POWERING THE UNIT & SELECTING OPERATING LANGUAGE AND REGIONAL PREFERENCES

- 1. Connect the power adapter to the rear panel of the meter.
- 2. Connect the power plug into the 24 V power socket.
- 3. Press the black ON/OFF power button.

At start up, the meter briefly displays the initialization screen.

4. The instrument launches into a Tutorial.

Tap \_\_\_\_\_ to automatically enter measurement screen.

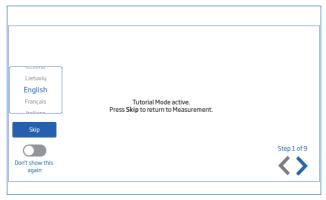

At first power on default language is English. A language window allows users to set operating language. To configure regional preferences, from Tutorial screen:

- Tap  $\equiv$  (direct **Menu** key) twice to access System Menu screen.
- Tap 
   (System Settings icon) and select System tab.
   Users can change the date, time, and region settings, as well as language preferences.

# 7.5. BASIC OPERATIONS

General operating modes are setup, measurement, logging, and data sharing.


- Tap  $\equiv$  (direct **Menu** key) to access:
  - 😬 User settings
  - System settings
  - O Measurement settings
  - Log recall files and file management

User can view a single sample or an interval log session, see Logging section for detailed description.

- Help for text and video support
- Tap 🔾 (direct **Home** key) to return to measurement.
- Tap 🗢 (Measurement Menu icon) to access sensor-related functions.

# 8. SYSTEM MENU ITEMS

Tap  $\equiv$  (direct **Menu** key) to access System Menu screen. Users must have previously logged in.

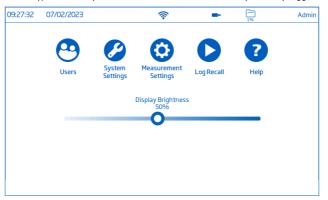

## System Menu capacitive icons

Symbol Name		Functionality		
•	Output         Users         Login and rights configuration & instrument accessibility			
Ø	System Settings	System configuration, connectivity, and printing items		
0	Measurement Settings	Measurement configuration		
0	Log Recall	Access logged measurement data		
?	Help	Access video-supported outline of main instrument functionalities		

## Brightness control bar



## 8.1. USERS

**Users** is the first item under the System Menu and enables multiple-users configuration, rights management, and account creation.



On first access, "Admin" is default user name and no password is required. Default options are updated from the Users menu.

Function	Administrator*	Standard User
Enable account creation	✓	_
Reset password	✓	_
Delete account	✓	_
Assign administrator rights	✓	_
View/Use Factory Settings reset	✓	_
Customize settings	✓	✓
Add FTP information	✓	✓
Change password	✓	✓
View and delete log files	✓	✓

\*Settings and configurations made by users with administrator rights can only be modified by users with identical rights.

#### Log in & Create New Account

- 1. Power the instrument. Wait for the initialization process to complete.
- 2. Tap to enter Measurement screen.

Note: Tap 🔹 (under 🔜) to disable Tutorial.

- 3. Tap  $\equiv$  followed by  $\mathfrak{S}$ .
- 4. Tap Edit Users to enter Account Management.

18:18:41 11/11/2022			(îr	<b>•</b> •	0%	HI6321
			User Settings			
Logou	t	Edit Users				

5. Tap  $\frown$  to enable Account Creation and Logins. Tap < to return.


<b>■</b> ■ 0%	HI6321
Enable Account Creation	
Enable Logins	

6. Tap Logout to enter Users screen.

"Admin" account is automatically created (default).

- 7. Tap the **plus** symbol avatar.
- 8. Input user name and tap 💌.

9. Enter password and tap 🛹. Reenter password to confirm.

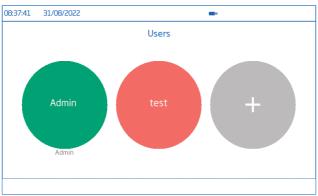

## Log Out & Switch User

- 1. Tap 😬 followed by Logout.
- 2. Tap user's account avatar.
- 3. Input password.

#### Adding & Removing Users (users with administrator rights only)

- 1. Tap  $\equiv$  followed by  $\mathfrak{S}$ .
- 2. Tap Edit Users to enter Account Management screen and start editing.

18:18:41 11/11/2022		((-	•	036	Admin
		User Settings			
Logout	Edit Users				
User Name			Admin		
Password			*****		
Icon Color					
Full Name			John Smith		
Info 1			Lab 01		

#### Configure user settings

Name, Password, assigned avatar color, Full Name, information fields, FTP-dedicated fields, Email Address

- To edit option, tap field and use the on-screen keypad to input information.
- Use the FTP dedicated fields and email address for file transfer of logged data.

#### Account Management

Users with administrator rights are able to:

- Enable Account Creation
- Enable Logins

Each power up requires user selection before instrument enters measurement mode.

- Enable Admin i.e. assign administrator rights to a standard user
- Reset Password
  - 1. Select user name from users list.
  - 2. Tap Reset Password.

The password is deleted and user can set a new password.

13:13:15	03/03/2022		<b></b>	<b>—</b> •	НІ6321
13.13.13	03/03/2022		<b>?</b>		0%
				Enable Accou	nt Creation
				En	able Logins 🔵
HI6321		Admin		Reset Password	Delete
user1		Admin		Reset Password	Delete


- Delete users
  - 1. Select user name and tap **Delete**. The instrument prompts for confirmation.

			1 1 M S	036	HI6321
HI6321 user1 user2		About to Delete	er? All associate	ed logs and	
	Confirm			Cancel	

## 8.2. SYSTEM SETTINGS

System Settings is the second item under the System Menu.

**Network**, **Connect** & **Print**, **System** tabs permit users to navigate system settings and operations, configure network connection and architecture, connectivity and printing services, change system settings, and view meter information.

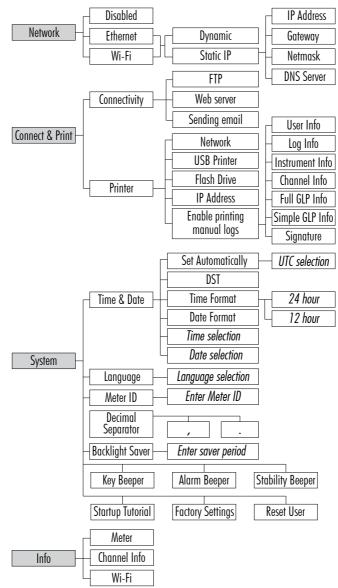

# Network

## Data sharing options: Ethernet, Wi-Fi, or Disabled

With connection established, IP assignment can be set as:

- Dynamic IP Address, Gateway, Netmask, DNS Server are auto assigned
- Static network details are entered manually

To input network information:

- 1. Tap IP Address field.
- 2. Input address and tap <
- Wi-Fi connectivity
- 1. Tap Wi-Fi.
- 2. Select the IP address type (Dynamic or Static).
- 3. Tap Press to select network.
- 4. Scan options and select preferred network. Enter password if / when prompted.
- 5. Tap < to confirm.

		System Setting	s		
Network	Disabled	Ethe	ernet	Wi-Fi	
onnect & Print	Dynamic	Static	Press to select	network	
System	IP Address				
Info	Gateway				
	Netmask	255.255.255.0			
	DNS Server	8.8.8			
	C C	,			

Note: With connection established, tap ♣ or <a>
 for to check IP address or verify connection status. When attempting to connect, <a>
 for is displayed <a>
 for sec. <

# **Connect & Print**

# **Options: Connectivity, Printer**

# Connectivity

Tap **connectivity** options:

- FTP access to meter: log file transfer to an FTP site and meter FTP server connection to client (log download)
- Meter web server: log file download to a web client
- Sending emails: log file transfer via email. A valid e-mail address is required for file sharing (see Users, Configure user settings).

18:47:35 11/11/2022		((;-	•	0%	HI6321
		System Settings			
Network	Connectivity				
Connect & Print	Allow F	TP access to meter			
System	Enable I	meter web server			
Info	Enable :	sending emails			
	Printer				
	Standard For	mat Receipt Fo	ormat		
Network		USB Prin	nter	<b>Flash Drive</b>	
	IP Address				

#### Printer

## Supported printer types: Standard, Thermal

- Select Standard Format to print on an international (ISO) or North American (ANSI) paper size standard.
- Select Receipt Format to configure information to be printed on a receipt format.

#### Print document options: Network, USB Printer, Flash Drive

- Select Network to connect a printer in the same network. Tap to enter IP address.
- Select USB Printer to connect a printer via USB-A port.
- Select Flash Drive to export log files directly to USB Flash Drive.
- Tap Enable printing manual logs to send files to configured printer.

Tap to enble for printing: User, Log, Instrument, and Channel information; Full GLP, Simple GLP, Signature.

**Note:** The connection to the printer (Flash drive) must be made prior to enabling the field selections. Ensure internet connection established between printer and internet.

18:49:43 11/1	1/2022	(îr	<b>—</b> ¤	0%	HI6321
		System Settings			
Network	Select the fie	lds to be printed			
Connect & Print		) User Info			
System		) Log Info			
		Instrument Info			
Info		Channel Info			
		Full GLP Info			
		Simple GLP Info			
		Signature			

# System

Options: Time, Date, Language, Meter ID, Decimal Separator, Backlight Saver, Beepers, Startup Tutorial, Factory Settings, Reset User

Note: Use the scroll bar to view or select from entire settings list.

# Time & Date

Tap **(**disable):

- Set Automatically (meter must be connected to the internet)
  - Direct selection from scrollable list of options
  - UTC options: from UTC 00:00 to UTC + 14:00 from UTC 00:00 to UTC -12:00



• DST (Daylight Savings Time) seasonal time change is used in some locations that advances clocks (typically by one hour) during warmer months.

09:57:14 03/03/2022		ŕ	<b>—</b> ¤	0%	HI6321
		System Settings			
Network	Time & Date	Set Automatically		UTC+01:00	
Connect & Print		DST	UTC	UTC+02:00	
System	J			UTC+03:00	
		Hour Minute Second		Time Format	
Info	Time	$ \begin{array}{c} 8 \\ 9 \\ 10 \end{array} : \begin{array}{c} 55 \\ 56 \\ 57 \end{array} : \begin{array}{c} 47 \\ 48 \\ 49 \end{array} $	AM PM	24 Hour 12 Hou	ır
	Date	Year Month Day 2021 2022 2022 2023 2023 2023 2023 2024 2 2 2 2 2 2 2 2 2 2 2 2 2		Date Format	
		Cestina			

# Time

• Hour, Minute, Second, time of day (AM or PM), time format (24 or 12 Hour) *Note:* Set Automatically must be disabled

Date: Year, Month, Day, choice of display formats ( DD-Mon-YYYY; YYYY-Mon-DD; DD/MM/YYYY; MM/DD/ YYYY; YYY/MM/DD; YYYY-MM-DD; Mon DD, YYYY )

Note: Set Automatically must be disabled

Language: select from list of supported options to change meter's interface language

#### Meter ID

- Name the meter with a discrete name, location, or number.
- Tap < to save.

Note: option available to users with administrator rights only.



Tap or corresponding tab to enable (disable) following settings:

Decimal Separator: comma or period depending on regional preference

Backlight Saver: enabled, 1 to 60 minutes (or disabled)

If the backlight turns off after the set period of time, tap screen to turn it back on.

Beeper: Key, Alarm, Stability

When enabled, an audible signal alerts users in the event of a wrong key press, an alarm condition, or the stability threshold being exceeded.

Startup Tutorial: if disabled (option enabled as default), the meter does not launch into Tutorial at power on.

12:01:05 20	23-02-16		-	• <u>1%</u>	Admin		
System Settings							
Network Connect & Print	Backlight Saver		10	minutes			
System		Key Beeper					
Info		Alarm Beeper					
		Stability Beeper					
		Startup Tutorial					
	Factory S	Settings	Reset User				

#### **Factory Settings**

Option restores system settings i.e. resolution for measured data, temperature unit, view mode, and alarm to original factory values. Restoring factory settings deletes all user information, logs, or configured measurement profiles. When option invoked, the instrument asks for confirmation.

Note: Option available to users with administrator rights only.

10:10:07	25/02/2022	(îr	<b>a</b> n <u>0%</u>	HI6321
Netwol	Re	store to Factory S	settings	
Systen	Are you su	re that you want to restore	to Factory Settings?	
	Confirm		Cancel	

#### **Reset User**

Option restores default settings for this user. All data (including profiles and log files) specific to this user will be permanently deleted, except for the username and password.

When option invoked, the instrument asks for confirmation.

#### Info

Read-only item displays information on meter, channel serial number, and Wi-Fi firmware version.

09:59:28 03	/03/2022	Ŕ	<b>.</b>	0%	HI6321
		System Settings			
Network Connect & Print	Meter:	Code: HI6321 Serial Number: 123456789LMN Firmware Version: 1.00 MAC Address: 70:1E:68:80:04:05			
System Info	Channel Info:	Type: EC Serial Number: 000214 Firmware Version: 1.00 Factory Calibration: 09/08/2022			
	Wi-Fi:	Firmware Version: 19.6.1			
	L				

# 8.3. MEASUREMENT SETTINGS

**Measurement Settings** is the third item under the System Menu and allows setting calibration, reading, temperature, view, alarms, logging and measurement profile options.



# **Measurement Setup Overview**

**Calibration, Reading, Temperature, View, Alarms, Logging, Profile** tabs help user navigate through all measurement operations. The following table presents an overview of possible functions.

## Calibration

	Candudidu	Desistivity	тос	Salinity	
	Conductivity	Resistivity	TDS	%	ppt, PSU
Last Calibration	Calibrate Clear	Based on conductivity or	Based on conductivity	Calibrate Clear	Based on conductivity
	ciour	salinity calibration	calibration	ciour	conductivity
Standard Entry Type		Automo	atic, Manual		
Calibration Reminder	Disabled Daily Periodic	-	_	Disabled Daily Periodic	-
Cell Constant	0.0500 to 200.0000	_	-	-	
Calibration Points	Single Point Multiple Points	_	_	_	

Reading

	Conductivity	Resistivity	TDS	Salinity			
	μS/cm	Ω∙cm	ppm	ppt			
Unit	mS/cm	kΩ∙cm	ppt	PSU			
Salinity Scale	AutoRanging	mΩ∙cm	AutoRanging	%			
		AutoRanging					
TDS Factor	-	-	0.40 to 1.00	-			
Stability Criteria		Accurate,	Medium, Fast				
Reading Mode		Direct					
Keuunig Moue		Direct/Autohold					
Probe Type		Pro	be code				

Temperature

	Conductivity	Resistivity	TDS	Salinity		
Temperature Source		Automatic, Manual				
Temperature Unit		°C, °F, K				
Temperature Compensation		Linear, Natural, Standard, Disabled				
Manual		-20.0 °C to 120.0 °C				
Reference Temperature	(whe	5.0 °C to 30.0 °C (when temperature compensation option is not disabled)				
Temperature Coefficient	(M	0.00 to 10.00 %/°C (with Linear temperature compensation option only)				
User Temperature Calibration	Calibrate Clear					

# View

	Construction	Destart in	TDC	Salinity	
	Conductivity	Resistivity	TDS	%	ppt, PSU
	Basic	Basic	Basic	Basic	Basic
	Simple GLP	Simple GLP	Simple GLP	Simple GLP	Graph
View Type	Full GLP	Graph	Graph	Full GLP	Table
	Graph	Table	Table	Graph	
	Table			Table	

# Alarms

		Conductivity	Resistivity	TDS	Salinity		
	High / Low	0.000 to	1.000 to	0.000 to	0.000 to 8.000E+1 ppt 0.00 to 42.00 PSU		
		1.100E+3 mS/cm	$1100E + 8 \Omega \bullet cm$	4.000E + 5 ppm	0.0 to 400.0%		
	High / Low	-20.0 to 120.0 °C					
	Temperature	-4.0 to 248.0 °F					
		253.2 to 393.2 K					

# Logging

	Conductivity	Resistivity	TDS	Salinity		
		Automatic Manual				
Logging Type						
	Autohold (Direct/Autohold mode only)					
Sampling Period		1, 2, 5, 10, 30 sec.				
(Automatic logging)		1, 2, 5, 10, 15, 30,	60,120, 150, 180	min.		
File Name		Create (Manua	l logging type only)			
Log Note		Us	er input			
Log Info 1 through to 4		Us	er input			
			None			
Sample ID	Increment					
	Manual					

# Profiles

	Conductivity	Resistivity	TDS	Salinity		
	Enable					
Profile Feature		Disable				
		Save As				
Current Profile		Save				
	Delete					
Load Profile		Configured profiles				

# 8.4. LOG RECALL

**Log Recall** is the fourth item under the System's Menu and allows data selection, viewing, sharing, and deletion. Logged data is retrieved by the user that has logged the measurement.



- Data is stored in parameter-specific .CSV files: conductivity, resistivity, TDS, salinity.
- Storage location is independent and organized in lots.
- A lot (file) can store 1 to 50 000 log records i.e. saved measurement data points.
- One user can store a minimum of 1 000 000 data points.

#### View

Data can be viewed, plotted (graph), or tabulated (complete with date, time, notes).

From the System Menu screen:

 Tap O (Log Recall). The instrument displays the Log Recall screen. Tap on corresponding table header element to order logged data by Name, Parameter, Start/Stop time stamp, #Samples.

Tap the up arrow to reverse order.


- 2. Tap to select .CSV file.
- 3. Tap View.

Note: Option available with .CSV file selected only.

04:21:34	11/11/2022	2	(	<b>a</b>	1%	HI6321
View	Select All	Deselect All	Log Reca	II	Delete	Share
	Name		Parameter		Start/Stop	#Samples
20221110_12	2913-ec_auto.c	:sv	Conductivity		9:13 10/11/2022 9:19 10/11/2022	7
20221110_12	2932-res_auto.	CSV	Resistivity		9:33 10/11/2022 9:46 10/11/2022	14
20221110_12	23021-tds_auto	.csv	TDS		0:21 10/11/2022 0:31 10/11/2022	11
20221110_12	23043-sal_auto	.csv	Salinity		0:43 10/11/2022 0:50 10/11/2022	8
20221110_12	23106-ec002.0	csv	Conductivity		1:06 10/11/2022 1:24 10/11/2022	10
20221110_12	23157-ec_Lab02	2_003.csv	Conductivity		1:57 10/11/2022 2:11 10/11/2022	9
				15:3	3:13 10/11/2022	

4. Tap 🗐 or 🔄 icon to have logged data displayed in tabulated form or plotted.

				(			1%	
0221110	)_122913-ec_au	ito.csv						
Index					EC abs		T[°C] ATC	
1	10/11/2022	15:29:13	0.000	µS/cm	0	µS/cm	25.0	"H"
Z	10/11/2022	15:29:14	0.000	µS/cm	0	µS/cm	25.0	"H"
З	10/11/2022	15:29:15	0.000	µS/cm	0	µS/cm	25.0	"H"
4	10/11/2022	15:29:16	0.000	µS/cm	0	µS/cm	25.0	"H"
5	10/11/2022	15:29:17	0.000	µS/cm	0	µS/cm	25.0	ОК
6	10/11/2022	15:29:18	0.000	µS/cm	0	µS/cm	25.0	ОК
7	10/11/2022	15:29:19	0.000	µS/cm	0	µS/cm	25.0	ОК

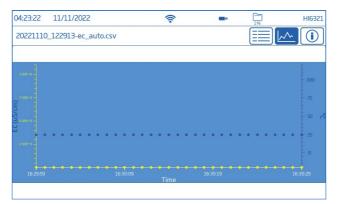

Tap 💿 icon and scroll through USER, LOG, CHANNEL, INSTRUMENT, GLP DATA information.

04:24:00 11/11/2022	(	<b>—</b> ¤	1%	HI6321
20221110_122913-ec_auto.csv				<u>~ ()</u>
LOG INFO				
Log Note:				
Log Info 1:				
Log Info 2:				
Log Info 3:				
Log Info 4:				
Log Type: Automatic				
Number Of Samples: 00000007				
Start Date and Time: 15:29:13 10/11/2020				
Sampling Period: 1 seconds				
Temperature Compensation: Linear				
Reference Temperature: 25.0°C				
Linear Compensation Coefficient:				
				_

# Select (Deselect) All

To export to a USB-A flash drive:

- 1. Tap (Log Recall) to access the log history.
- Tap Select All button to select entire log history. With all files selected, tap Delete to empty the log or tap Share to transfer data.
- 3. Tap **Deselect All** to clear selection.

#### Delete

- 1. Tap (Log Recall) to access the log history.
- 2. Tap to select required .CSV file (one file) or tap Select All.
- 3. Tap Delete.
- The instrument prompts for confirmation. Deleted files can not be recovered and Log Recall screen is displayed blank.

04:21:34	11/11/2022	(îr	<b>••</b> 1%	HI6321
View	Select All Deselect All	Log Recall	Delete	Share
	Del	ete Selected Log Fil	es?	amples
20221110_3				7
20221110_	Are you sure you	want to delete the selected log be undone.	g files? This cannot	14
20221110_				11
20221110_:				8
20221110_				10
20221110_	Confirm		Cancel	9
			15:33:13 10/11/2022	

# Share

# Options: USB-A, FTP, Email, Print, Web server USB-A and USB-C

Plug the USB flash drive into the USB port located on the back.

- 1. Tap (Log Recall) to access the log history.
- 2. Select file (files) for transfer or use Select All option.
- 3. Tap Share. Pop-up flyout is displayed.
- 4. Tap to select USB-A or USB-C.
  - $\overline{\mathbf{x}}$  is displayed during data transfer.
- 5. Transfer completion is confirmed and the instrument returns to Log Recall screen.

# FTP

HI6321 can act as an FTP server (host) or client. Meter has to be connected to the internet and Allow FTP access to meter enabled. See System Settings section, Connect & Print tab.

- Use meter's IP address and password to connected and view logged files.
- Enter in the FTP dedicated fields own server information to export logged files to the FTP server.
- Configure FTP server info in the User menu (😁) to use the meter as an FTP client and upload files to an FTP server.

# Connect via FTP to meter server:

- 1. On preferred FTP software, type the meter's IP address in the dedicated Host field.
- 2. Enter the username and password of the user currently logged in.
- 3. Connect to view the files logged on the meter.



Connect the meter to an FTP server and share logs:

- 1. In Users menu, type selected server's IP address, username, and password.
- 2. From System Menu, tap 오 . Log Recall screen opens up.
- 3. Select file(s) for transfer.
- 4. Tap Share. Pop-up flyout is displayed.
- 5. Tap to select FTP. The files are being transferred in the root folder of the server.

# FTP server installation and configuration

- PC running Windows10 or later
- Password protected Windows account
- FTP server must be allowed through the Windows Firewall

Installation

- 1. Navigate to Start > Control Panel > Administrative tools > Server Manager.
- 2. Go to Roles and expand Web Server.
- 3. Right click on Web Server and then click on Add Role Services.
- 4. Go to Role Services and check FTP Server.
- 5. Ensure IIS Manager (Internet Information Services) is checked under Management Tools.
- 6. Click Next followed by Install.
- 7. Wait for installation to complete.

Configuration (PC must be running Windows10 or later)

- 1. Navigate to Start > Control Panel > Administrative tools > IIS Manager (Internet Information Services).
- 2. Double click to expand the IIS Manager console.
- 3. Right click on Sites, on the Connection pane.
- 4. Click on Add FTP Site, to select. Type the FTP server name and the path to be used for file transfer *Note:* Select Make New Folder to create a designated folder to store FTP files.

- 5. Click Next.
- 6. In the Binding and SSL Settings window keep all default settings but change the SSL option to No SSL.
- 7. Click Next.
- 8. When prompted to authenticate and authorize information, select Basic and Specified users.
- 9. Type local account name to gain access to the server.
- 10. Check both **Read** and **Write** options.
- 11. Click **Finish**.

## Email

Meter has to be connected to the internet and **Enable sending emails** enabled (see System Settings, Connect & Print section).

Tap 🕙 to access the Users menu and input email address.

18:18:41 11/11/2022	🛜 📼 📴 Hi632:
	User Settings
Info 3	info text 3
Info 4	info text
Email Address	user.HI6321@email.com *required for file sharing
FTP IP Address	10.168.19.10
FTP User Name	ftp-user-name
FTP Password	*****

- 1. From System Menu, tap 오 to access Log Recall.
- 2. Select file(s) for transfer or use Select All option.
- 3. Tap Share. Pop-up flyout is displayed.
- 4. Select Email.
- 5. Transfer completion is confirmed and the instrument returns to Log Recall screen.

View	Select All Deselect A	Log Recall	Delete	Share
	Name	Parameter	Start/Sto <sup>Share T</sup>	io: les
	2012	6 . I. I. I.	15:29:13 10/11	USB-A
20221110 12	2913-ec_auto.csv	Conductivity	15:29:19 10/11	FTP
00001110 10	2022	Desistivity	15:29:33 10/11	
20221110_122932-res_auto.csv		Resistivity	15:29:46 10/11	Email
	2021 14	TDS	15:30:21 10/11	Print
20221110_123021-tds_auto.csv		TDS	15:30:31 10/11	
	2042	<b>C</b> (1) (1)	15:30:43 10/11	Cancel
20221110_12	3043-sal_auto.csv	Salinity	15:30:50 10/11/2022	
00001110 10	2105 002	Construction	15:31:06 10/11/2022	10
0221110_12	3106-ec002.csv	Conductivity	15:31:24 10/11/2022	10
	2157	Constant No.	15:31:57 10/11/2022	0
20221110_123157-ec_Lab02_003.csv		Conductivity	15:32:11 10/11/2022	9
			15:33:13 10/11/2022	

#### Print

- Connect either a printer (Network or USB) or plug-in a USB Flash Drive (see System Settings, Connect & Print section).
- Tap Print and follow on-screen instructions.

#### Web server

Any browser can be used to access the web server and download log files. Meter has to be connected to the internet and **Enable meter web server** enabled (see System Settings, Connect & Print section).

Note: Both the meter and the device the browser is accessed from have to be connected at the same network.

1. Tap 🛜 for IP address and type address in the browser.



2. Enter the username and password of current user to gain access to logs and tags. Click on file to download to the PC.

Log	in		System Information			
Username	Password		Category	System Information		
HI6321 Login	••••	1~	Meter:	Code: HI6321 Serial Number: 123456789 Firmware Version: 0.1.221209 RD1 MAC Address: 70:1E:68:80:00:01		
			Channel Info: Type: EC Serial Number: ( Firmware Versio Factory Calibrati		n: 0.2.7	
			Wi-Fi:			
			Wi-Fi:			
HANNA Instruments GS	Parameter	Number of Entries		Ston Time	1	
<b>gs</b> ille Name	Parameter Conductivity	Number of Entries	Wi-Fi: Start Time 20:32:17 12/12/22	Stop Time 22 20:32:24 12/12/2022	Download	Vie
GS ile Name 10221212_193217-ec_auto.csv	Conductivity		Start Time	20:32:24 12/12/2022	Download	
<b>GS</b> ile Name 10221212_193217-ec_auto.csv 10221212_193249-ec_Lab01_002/	Conductivity csv Conductivity	8	Start Time 20:32:17 12/12/20	20:32:24 12/12/2022           20:33:00 12/12/2022		Vie Vie
US IIIe Name 0221212_193217-ec_auto.cev 0221212_193249-ec_Labo1_002. 0221212_193321-ec_Labo1_003.4	Conductivity csv Conductivity	8	Start Time 20:32:17 12/12/20 20:32:51 12/12/20	20:32:24 12/12/2022           20:33:00 12/12/2022           20:33:38 12/12/2022	Download	Vie
	Conductivity Conductivity Conductivity Conductivity	8 10 9	Start Time 20:32:17 12/12/20 20:32:51 12/12/20 20:33:25 12/12/20	20:32:24 12/12/2022           20:33:24 12/12/2022           20:33:00 12/12/2022           20:33:38 12/12/2022           20:33:44 0 12/12/2022	Download Download	Vie

# PC Connection

The logged data can be transferred from the meter to a PC.

- Use the USB-C cable to connect the meter to the PC.
- The meter will appear as a flash drive on the computer.
- Save files to the PC. All logs will be listed as .CSV files.

The .CSV files may be opened with any text editor or spreadsheet application.

All features of the spreadsheet program can be used to analyze and graph the data.

## 8.5. HELP

Help is the fifth item under the System Menu.

• Tap 7 to access support and navigate through an overview of system's main functionalities.

09:27:32 07/02/2023		1%	Admin
Users		Stem Settings	
		Tutorial	
1. HI6321 First Look	$\sim$	1.2. Meter Layout	
1.1. Features		Front view	
1.2. Meter layout		Stability Measurement Start Measure	
1.3. Specifications		Indicator Profile Logging Setti	ngs
1.4. User Interface 1.5. Setting up the device	$\mathbf{\bullet}$	ANNA	
		Biotan_ii	
2. Getting Started	>	1416 <sup>µS/cm</sup>	
3. General Operations	>	25.0 %	
4. Logging	>		
5. Troubleshooting guide	>	< ○ ≡	
6. Accessories and Warranty	>		
		Back Home Menu	

- Tap to play (stop) video-supported segments:
  - ▶ 1.5. Setting up the device
  - ▶ 3.1. Calibration
  - ▶ 3.2. Reading measurements
- Tap  $\equiv$  once to increase video speed.
  - ▶ The meter supports, with each tap, three playback speeds: normal (x1), medium (x2), and fast (x4).

# 9. MEASUREMENT & PROBE SETUP MENU

From the Measurement screen, tap  $\heartsuit$  to access system and measurement configuration tabs. Alternatively, tap  $\equiv$  (**Menu** key) then O.



## 9.1. CALIBRATION

**Options: Conductivity & Salinity** 

### Last Calibration, Standard Entry Type, Calibration Reminder, Cell Constant, Calibration Points

Last Calibration (Conductivity and Salinity): Calibrate or Clear a previous calibration

#### Standard Entry Type

- Automatic: meter selects the closest calibration standard value to that of the sample being measured
- Manual: user selects the calibration standard to be used for calibration

Calibration Reminder (Conductivity and Salinity)

- Daily-calibration interval (hours and minutes)
- Periodic-calibration interval (days, hours, minutes)

"Calibrate probe" message is displayed if calibration is not done within set interval.

06:15:46 1	1/11/2022	ŝ	<b>•</b> •	HI6321						
	Measurement Settings									
Calibration	Last Calibration	Calibrate	Clear							
Reading Temperature	Standard Entry Type	Automatic	Manual	]						
View	Calibration Reminder	Disabled	Daily	Periodic						
Alarms			Hour Minute	Days Hours Minutes						
Logging	Cell Constant	1.0000	/cm							
Profiles	Calibration Points	Single Point	Multiple Points							

Cell Constant: allows users to adjust cell measurement to compensate for the measured properties.

- Tap dedicated field and enter value.
- Tap 🕶 to save.

Calibration Points: allows users to select a Single Point or Multiple Points calibration.

10:51:22	04/04/2022	2		((:-			D.	0%		HI6321	
	Measurement Settings										
Calibration	Last Cal	bration	Calibrate Clear								
Reading	Calibri	Enter c	ell cor	nstant value		1	z	з			
Temperatur View	e Calibra Remir	Range: 0.0500 ↔ 200.000 /cm				4	5	6			
Alarms			1.00			7	8	9	◀	es	
Logging	Cell Co	Cancel Save				0		-	Exp		
Profiles	Calibrati	on Points	Si	ingle Point	М	ultiple P	oints				

#### 9.2. READING

Options: Parameter, Units/Salinity Scale, TDS Factor, Stability Criteria, Reading Mode, Probe Type

16:51:46 20	023-02-23	ŕ		<b>—</b> •	2%	HI6321
		Measurement	t Settings			
Reading	Parameter	meter Conductivity Resistivity TDS			Salinity	
Temperature	Unit	ppm ppt			AutoRanging	
Alarms	View Alarms					
Logging	Stability Criteria	Accurate	te Medium		Fast	
Profiles	Reading Mode	Direct	Direct/A	utohold		
	Probe Type	HI7631233				

#### Parameter

Options: Conductivity, Resistivity, TDS, Salinity

Tap to select measurement configuration.

#### Units & Salinity Scale

- Conductivity Options: µS/cm, mS/cm, AutoRanging i.e meter automatically selects the unit to optimize the measurement.
- Resistivity Options: Ω·cm, kΩ·cm, MΩ·cm, AutoRanging Resistivity is calculated from conductivity measurement.

- TDS Options: ppm, ppt, AutoRanging i.e. meter automatically selects the unit to optimize the measurement.
- Salinity scale: ppt (Natural Sea Water 1966), PSU (Practical Scale 1978), %(Percent Scale) Salinity value is determined based on selected scale.

## TDS Factor

TDS factor is a conversion factor used to change an EC measurement to a ppm measurement. TDS factor may be set between 0.40 to 1.00.

- Tap dedicated field and enter the value.
- Tap < to save.



# Stability Criteria

# Options: Accurate, Medium, Fast

Allows users to set the stability criterion for selected parameter based upon the stability setting selected. While the measurement is changing, the stability indicator is shown partially (\_\_\_\_\_\_); when the criteria is reached the indicator is displayed as a full green bar (\_\_\_\_\_\_).

- Accurate: for applications where high accuracy is required Measurement is recognized as stable using more critical criteria evaluating measurement fluctuations.
- Medium: for applications where average accuracy is accepted Measurement is recognized as stable using less critical criteria evaluating measurement fluctuations. The measurement may still change after registering stable.
- Fast: for applications when stability-evaluation methods are not so critical

# **Reading Mode**

# Options: Direct, Direct/Autohold

- Direct: as measurement changes, measurement stability is continuously evaluated. Unstable (blinking) or Stable indicator are displayed above the progress bar.
- Direct/Autohold: 1 (lock icon) is displayed between the log and Measurement Menu icons.
  - ▶ Tap to initiate a measurement.

- ▶ Configured parameter icon is displayed i.e. 🕫 📼 📼 🖘.
- \_\_\_\_\_ is displayed blinking until the measurement is stable.
- ▶ When stable, the indicator stops blinking and measurement freezes at current value.
- ▶ Tap parameter icon to return to direct measurement.



#### 9.3. TEMPERATURE

Options: Temperature Source, Temperature Unit, Temperature Compensation, Manual, Reference Temperature, Temperature Coefficient, User Temperature Calibration

06:21:28 11	1/11/2022	🔷 🗣 🗎			HI6
		Measureme	ent Settings		
Calibration Reading	Temperature Source	Auto	matic	Mar	nual
Temperature	Temperature Unit	°C	c	F	К
View	Temperature Compensation	Linear	Natural	Standard	Disabled
Alarms Logging	Manual	25.0	°C		
Profiles	Reference Temperature	25.0	°C		

#### Temperature Source: Automatic, Manual

Users can select between physical temperature input source (Automatic) and entering sample temperature value manually (Manual).

- Automatic: built-in temperature sensor adjusts the measured conductivity to a reference temperature. "ATC" is displayed indicating temperature compensation status.
- Manual: sample temperature is entered manually and displayed "MTC" indicates temperature compensation status.

### Temperature Unit: Celsius, Fahrenheit, Kelvin degrees

### Temperature Compensation (Scale): Linear, Natural, Standard

- Linear: used when it is assumed that the temperature coefficient of variation has the same value for all measurement temperatures.
- Natural: for natural ground, well, or surface water and covers 60 to 1000  $\mu$ S/ cm (0 to 35 °C) range.
- Standard: used for resistivity measurements (high-purity water).

### Manual

Manually input sample's temperature when Temperature Source is configured as Manual.

- Tap dedicated field and enter the value.
- Tap < to save.

# Reference Temperature: 5.0 to 30.0 °C

Value used for temperature-compensated conductivity. All conductivity measurements will be referenced to the conductivity of a sample at this temperature.

- Tap dedicated field and enter the value.
- Tap < to save.

# Temperature Coefficient: 0.00 to 10.00 %/°C (Linear Scale only)

Temperature coefficient is a function of the solution being measured. When the actual temperature coefficient of the sample is known:

- Tap dedicated field and enter the value.
- Tap < to save.

## User Temperature Calibration: Calibrate or Clear

To perform a new calibration:

- 1. Tap Calibrate.
- 2. Place the probe and a reference thermometer into a stirred container of water. Allow reading to stabilize.
- 3. If the displayed value is different than the reference thermometer reading, tap **Calibrated Temperature** and use the on-screen keypad to enter the value.

Note: Users need a separate reference thermometer.

- 06:21:40 11/11/2022 <u></u> -HI6321 Measurement Settings Temperature Natural Standard Disabled Compensation Manual 25.0 °C Temperature Reference 25.0 °C Temperature Temperature Coefficient 1.90 %/°C Last Calibration User Temperature Calibrate Clear 11/11/2022 Calibration 22:39:35 11/11/2022 <u></u> HI6321 Calibrate temperature Measured Temperature: °C 25.0 °C **Calibrated Temperature:** Check if the displayed value is different than the reference thermometer reading. Touch the Calibrated Temperature value to edit value to reference temperature. Press 'Save' to complete calibration. Cancel Save
- 4. Tap Save to confirm and save the calibration data.

#### 9.4. VIEW

Conductivity, Salinity %: Basic, Simple GLP, Full GLP, Graph, Table Resistivity, TDS, Salinity ppt & PSU: Basic, Simple GLP, Graph, Table

• Select preferred display configuration from View Type window.

10:49:30 03/03/2022	4	(h.	<b>•</b> •	0%	HI6321		
Measurement Settings							
Calibration Reading	Basic Simple GLP						
Temperature View Type	Full GLP						
	Graph						
View	Table	J					
Alarms							
Logging							
Profiles							

#### Basic

Screen displays the measured value, measurement unit, stability and temperature compensation status/source.



### Simple GLP

Simple GLP screen displays last calibration date and time, offset value and cell constant (as well as data displayed in Basic screen).

Note: If no calibration was made, "Not Calibrated" is displayed.



### Full GLP

Full GLP screen displays probe symbol, calibration temperature, calibration information (as well as data displayed in Simple GLP screen).



#### Graph

When Graph is selected the measured value is plotted as a graph.

To zoom in on a graph:

- 1. Select Time or parameter axis.
- 2. With axis selected, pinch (drag) on the display.

1416 <sup>µS/cm</sup>	25.0°C	. 🕘 🚽	<b>P</b>
1410	LJ.OATC		<b>1</b>
			- 100
			- 75
			- - 50 -
⊇ soores_  			
			- 25
		TRACKING	

## Table

When Table is selected, the measured values are displayed tabulated (complete with date, time, and notes made during logging). The newest data is displayed at the top of the table.

14:01:38 31/08	3/2022		<b></b>	<b>••</b> 0%	HI6321
Stable 1416	µS/cm	25.0°C	rc	•	Ф
EC (µS/cm)	T(°C)	Time	Date	Notes	
1416	25.0	14:01:37	31/08/2022		
1416	25.0	14:01:36	31/08/2022		
1416	25.0	14:01:35	31/08/2022		
1416	25.0	14:01:34	31/08/2022		
1416	25.0	14:01:33	31/08/2022		
1416	25.0	14:01:32	31/08/2022		
1416	25.0	14:01:31	31/08/2022		
1416	25.0	14:01:30	31/08/2022		
1416	25.0	14:01:29	31/08/2022		
1416	25.0	14:01:28	31/08/2022		
1416	25.0	14:01:27	31/08/2022		

## 9.5. ALARMS

## Options: High /Low configured parameter value, High/Low Temperature

Users can set the threshold limits for the measured parameters. When the measurement exceeds set high-threshold value or drops below set low-threshold value, the alarm is triggered and an alarm message is displayed. If the Alarm Beeper is enabled (See: System Menu Items / System Settings / Alarm Beeper), an audible beep will be heard.

Note: High-parameter values cannot be lower than low-parameter ones.

To set an alarm limit:

- 1. Tap 🔵 to enable low or high parameter threshold option.
- 2. Use the on-screen keypad to enter the value.
- 3. Tap Enter to confirm (Cancel to exit and return to measurement settings).

12:45:43 03	12:45:43 03/04/2022 🛜 📼 🛱									
	Measurement Settings									
Calibration	High Conductivity		0.000	μS/cm						
Reading	Low Conductivity		0.000	μS/cm						
Temperature	High Temperature		50.0	_ ℃						
View	Low Temperature		10.0	°C						
Alarms										
Logging										
Profiles										

🔯 (alarm icon) is displayed on the measurement screen when an alarm is active.



#### 9.6. LOGGING

Options: Logging Type, Sampling Period, File Name, Log Note, Log Info, Sample ID

12:38:29	03/04/2022	ŝ	<b>—</b> ¤	0%	HI6321
	_	Measurement S	ettings		
Calibration	Logging Type	Automatic	Manual	Autohold	
Reading Temperature	Sampling Period	180 min 1 sec 2 sec			
View	File Name	Create	UnnamedLog		
Alarms	Log Note	Lab 112			
Logging	Log Info 1	test			
Profiles	Log Info 2				

### Logging Type

- Automatic: data is logged automatically at predefined time intervals (i.e. Sampling Period).
  - ▶ A file name is automatically generated, complete with year/month/day, and logging time.
  - Log files are identified by parameter:
    - EC: 20220329\_085101-ec\_auto
    - Resistivity: 20220309\_095704-res\_auto
    - Salinity: 20220331 101507-**sal** auto
    - TDS: 20220331\_101458-tds\_auto
- Manual: data is logged every time 🐠 is tapped.
  - Tap Create to rename file as preferred e.g. 20220329\_085101-ec\_Test\_001 or 20220329 085101-tds 002.
- Autohold: Direct/Autohold reading mode must be configured.

**Note:** Manual and Autohold records are stored in the same log file, i.e. data logged on different days is stored in the same lot. Automatic records are stored separately. Data logged with Autohold option selected, is identified by the "H" in the Notes column.

## Sampling Period

Option available only with **Automatic** logging type selected. Time-interval options are selected from scrollable list.

# File Name

Available only with Manual and Autohold logging type selected.

To create a file name, from Logging screen:

- 1. Tap Create.
- 2. Use the on-screen keypad to enter file name (maximum 13 characters).
- 3. Tap Enter to confirm.

# Log Note & Log Info

Notes on measured data are saved together with logged data.

# Sample ID

Manual and autohold samples can be labeled with a numerical ID (increments with each new sample logged), a text label, or a text label with numerical ID.

With Increment selected:

- 1. Tap Sample ID Prefix.
- 2. Use the on-screen keypad to enter a text prefix, of maximum 15 characters, to the sample name.
- 3. Tap < to confirm.
- 4. Select ID number from scrollable list.

# 9.7. PROFILES

Option: Profile Feature, Current Profile (configure), Load Profile (previously configured)

• Tap 💿 to enable (disable) Profile Feature option.

# **Current Profile**

A profile is a sensor setup complete with required measurement and temperature unit, display preference, and alarm threshold options. Once configured and saved the profile can be loaded for applications that require similar configurations.

To save a profile, having previously configured all other application-specific options, from the Measurement screen:

- 1. Tap 🗘.
- 2. Tap Profiles tab.
- 3. Tap Save As and use the keypad to enter profile name.
- 4. Tap < to confirm.

Once saved, profile name is added to the Load Profile list.

# Load Profile

To select a configured profile:

- 1. Tap to select from the Load Profile list.
- 2. Profile name is automatically entered in the Current Profile field.

3. Start measuring.

		0%	HI632
Measurement Settings			
Chem-Lab-02 (Modified)			
Save As S	Save	Delete	
Chem-Lab-02			
EC_daily			
	Chem-Lab-02 (Modified) Save As Chem-Lab-02	Chem-Lab-02 (Modified) Save As Save Chem-Lab-02	Chem-Lab-02 (Modified) Save As Save Delete Chem-Lab-02

Note: Profile setting modifications are indicated by (\*) in measurement screen.

# **10. CALIBRATION**

# **10.1. CONDUCTIVITY CALIBRATION**

HI6321 meter allows two types of calibration procedures:

- Standard up to four-point conductivity calibration with standards for cell factor determination:
  - ▶ 0  $\mu$ S/cm for offset
  - ▶ 84 µS/cm, 1413 µS/cm, 5.0 mS/cm,12.88 mS/cm for the 0.1/cm cell
  - ▶ 80.0 mS/cm, 111.8 mS/cm additional standards for the 1.0/cm cell
- Standard single point salinity calibration in 100% salinity standard, with the parameter reading set to Salinity and Salinity Scale set to %.

Note: Conductivity parameter must be calibrated first.

### Preparation

- Remove plastic bung prior to calibration.
- Clean the probe in distilled water, shake off water droplets, and allow to dry prior to calibration.
- Use a calibration standard with a value that is close to that of the sample.
- Inspect the probe for debris or blockages.
- Ensure the vent holes are completely submerged.
- Tap the probe to remove any air bubbles that may be trapped inside the sleeve.
- To minimize cross-contamination, when a two-point calibration is required, use two beakers: one for rinsing the probe and the other for calibration.
- For measurements across a temperature gradient (when water temperature is drastically different from the standards), allow the probe to reach thermal equilibrium before conducting calibrations or making measurements.

### 10.2. PROCEDURE

#### Automatic Calibration

With option selected, the HI6321 automatically selects the closest standard to the value being measured.

12:18:32	30/01/202	3		()	>	-		0%	HI6321
Stab	le 1	413 μ	ıS/cm	MTC	25.0	°C	Cell K.:	0.9942 /cm	
Calibrate Conductivity mS/cm Confirm standard Touch 'Confirm Standard' to update the calibration.								I	
Clea	ar Calibration		Cancel		Save			Confirm Standar	ď
	mS/cm 5.000 2.000 mS-> 20.0	J Lms							
	0.9942/cm 25.0 °C 30/01/202: 12:15:36								

With the probe connected to the meter:

- 1. From Measurement screen, tap 🗢 (Measurement Menu icon).
- 2. Tap Temperature tab.
- 3. Configure Temperature Source.
- 4. Tap Calibration tab.
- 5. Configure Standard Entry Type as Automatic.
- 6. Select number of calibration points.
- 7. Tap Calibrate to start the process.
- 8. Tap **Clear Calibration** to start a fresh calibration and clear all previously used standards. If using a new conductivity probe, it is recommended to clear all standards and start new.
- Suspend the probe in the air. Allow for the reading to stabilize. The standard value is automatically recognized.
- 10. After the reading has stabilized, tap **Confirm Standard**. The calibration point is added to the tray.


### Single Point calibration

Tap Save to have first calibration point accepted and saved. Meter returns to Channel Settings.

## **Multiple Points calibration**

Use two beakers, one for rinse and one for measurement.

- 11. Raise and lower the probe in rinse beaker of standard. Discard rinse standard.
- 12. Immerse the sensor in standard. The standard value is automatically recognized.
- 13. Allow reading to stabilize then tap Confirm Standard.
- 14. Repeat steps 11 through to 13 with up to two more standards.
- 15. Tap Save to accept and save (in tray) the calibration point and return to calibration settings.
- 16. Tap  $\bigcirc$  or < to return to measurement.

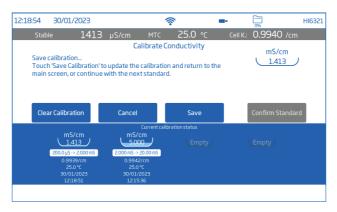

### **Manual Calibration**

With option selected, users have to option to edit calibration point values.

- 1. Follow Automatic Calibration procedure up to step 9 but configure Standard Entry Type (Step 5) as Manual.
- 2. Suspend the probe in the air. Allow for the reading to stabilize.
- 3. Tap Edit Standards. Input value and tap Save.
- 4. After the reading has stabilized, tap Confirm Standard.
- 5. Raise and lower the probe in next rinse standard. Discard.
- 6. Immerse the sensor in calibration standard. Tap Edit Standards.
- 7. Input value and tap Save.
- 8. After the reading has stabilized, tap Confirm Standard.
- 9. Continue with up to two more standards.



### **10.3. SALINITY CALIBRATION**

Salinity calibration can be performed with the parameter reading set to Salinity and Salinity Scale set to %.

- 1. From Measurement screen, tap 🗢 (Measurement Menu icon).
- 2. Tap Reading tab.
- 3. Tap to configure reading Parameter to Salinity.

- 4. Set Salinity Scale to %.
- 5. Tap Calibration tab.
- 6. Tap **Calibrate** to start the process.
- 7. Tap **Clear Calibration** to start a fresh calibration. If using a new conductivity probe, it is recommended to clear and start new.
- 8. Raise and lower conductivity cell in rinse beaker of salinity standard then discard.
- 9. Immerse the sensor in salinity standard. Ensure the vent holes are completely submerged. The meter automatically recognizes the standard.
- 10. Allow the reading to stabilize then tap **Confirm Standard**. The standard is recognized and added to the tray.
- 11. Tap Save to have calibration accepted and saved, and return to calibration setting screen.

06:36:58	15/11/2	2022		1	(h.	•		) нюз21
Stab		100.0			25.0			
Calibrate								
Maxin	Calibration complete Maximum number of calibration points reached. Touch 'Save Calibration' to update the calibration and return to the main screen.							
Clea	ar Calibrati	on	Cancel		Sav	e	Cor	nfirm Standard
			Current M	10D1 EC (	alibratior	n status		
	Lat Calibration Date 15/11/2022 00:06:19							

# **11. MEASUREMENT**

## Options: Direct, Direct/Autohold

- Direct: sample measurements are displayed continuously.
- Direct/Autohold: reading is displayed when measurement stability is reached. A measurement that has not reached equilibrium will not be used. After stability criteria is reached, the meter enters Direct/Autohold mode.

## When to measure conductivity instead of resistivity

Resistivity is the reciprocal of conductivity and their scales emphasize different areas of the measurement range.

- Resistivity (set as Reading parameter) is commonly used in ultrapure water
- Conductivity (set as Reading parameter) is suitable for measuring larger amounts of contaminants. Users can subsequently change Reading parameter to Resistivity to measure in resistivity units (MΩ • cm).

Recommended temperature compensation setting for these type of measurements is **Standard** as it utilizes the correct compensation algorithm.

## 11.1. MEASUREMENT TIPS

- 1. Connect the probe to the meter. Only work with recently calibrated probes.
- Use the HI764060 electrode holder for easy transfer in and out of containers during calibration and sample measurement.
- Ensure plastic bung is removed prior to taking measurements. To limit sample contamination, pour 2 beakers of calibration standards. Use one beaker to rinse the sensor, and another one for measurement.
- 4. Ensure the vent holes are completely submerged.
- 5. Tap the probe to remove any air bubbles that may be trapped inside the sleeve.

Note: Use the same size beaker and immersion depth for samples and calibration standards.

- 6. If measuring across a temperature gradient, allow the sensor to reach temperature equilibrium. If using manual temperature compensation, input the sample temperature.
- 7. Once the reading indicates Stable, record measurement data.

# 11.2. DIRECT READINGS

- From Measurement screen, tap 🌣 (Measurement Menu).
- Tap Reading tab and select measured parameter.
- Tap **Direct** to select direct Reading Mode.
- Place the probe into the sample to be measured. Ensure the vent holes are completely submerged. Allow time for the reading to stabilize. (Unstable status is indicated on the screen until measurement is stabilized).
- The measured value is displayed on the LCD, together with GLP information and display preferences.



# 11.3. DIRECT / AUTOHOLD READINGS

- From Measurement screen, tap 🌣 (Measurement Menu).
- Tap Reading tab and select measured parameter.
- Tap **Direct/Autohold** to select direct/autohold Reading Mode.
- Place the probe into the sample to be tested. Ensure the vent holes are completely submerged.

- Measured value is displayed on the LCD.
  - is displayed blinking until configured stability criteria is reached.

When measurement is stable, the measured value is kept on display and Automate stops blinking.

• To return to direct Reading Mode, tap icon on display ( @/@/@/).



# 12. LOGGING

Tap 🚞 to check available storage.

10:07:11	03/03/2022	()	e	HI632
Storage	Used: 0.07 MB / 25	5 84 MR		Close
Storage	03eu. 0.07 HD7 25	5.64140		Close

Three logging types are available: Automatic, Manual, and Autohold.

12:38:29 03	3/04/2022	<u></u>	<b>—</b> ¤	0%	HI6321
	Measurement Settings				
Calibration	Logging Type	Automatic	Manual	Autohold	
Reading Temperature	Sampling Period	180 min 1 sec 2 sec			
View	File Name	Create	UnnamedLog		
Alarms	Log Note	Lab 112			
Logging	Log Info 1	test			
Profiles	Log Info 2				

#### Automatic logging

- Readings are logged (
   ) at predefined period intervals. Sampling interval options range from 1 second to 180 minutes.
- Records are continuously added to it until the session stops.

- For each automatic logging session, a new lot is created.
- A complete set of GLP information including date, time, mode selection, temperature reading, and calibration information is stored with each log. User and log information e.g. company, address, sample details may also be included on the .CSV file.

## Manual logging

- Readings are logged each time 🜑 is tapped.
- All manual readings are stored in a single log (i.e. records made on different days share the same log).

# Autohold logging

- Readings are logged each time 🔍 is tapped and configured stability criteria is reached.
- All Autohold readings are stored in a single log (i.e. records made on different days are logged in the same lot).

Note: Stability criteria can be set to Fast, Medium, or Accurate for all logging types.

# 12.1. AUTOMATIC LOGGING

- 1. From Measurement screen, tap 🗢 (Measurement Menu).
- 2. Tap Reading tab and tap to select stability criteria (Accurate, Medium, or Fast).
- 3. Tap Logging tab and select Automatic logging type.
- 4. Scroll to select Sampling Period.
- 5. Tap 🔾 (Home key) to reenter Measurement screen.
- 6. From Measurement screen, tap 🗢 symbol to start/stop logging.
- 7. During active logging, users can track and log notes. To add a new log note:
  - Tap 📃 .
  - Use on-screen keypad to enter text.
  - Tap < to log entered note. Log Recall view, and .CSV log indicate the added log notes.

# 12.2. MANUAL LOGGING

- 1. From Measurement screen, tap 🗢 (Measurement Menu).
- 2. Tap Reading tab then select stability criteria (Accurate, Medium, or Fast).

**Note:** Manual logging doesn't use this criteria for logging. Setting this is only for the user to note the stability of the reading.

- 3. Tap Logging tab and select Manual logging type.
- 4. See option to create file name:
  - Tap Create.
  - Use on-screen keypad to enter file name.

• Tap < to save new name.

12:40:20 12	1/11/2022	(îr	<b>•</b> •	0%	HI6321
		Measurement	Settings		
Calibration	Logging Type	Automatic	Manual	Autohold	
Reading Temperature	Sampling Period	180 min 1 sec 2 sec			
View	File Name	Create	UnnamedLog		
Alarms	Log Note	Lab 112			- 1
Logging	Log Info 1	test			
Profiles	Log Info 2				

- 5. See option to define sample ID (name and prefix).
  - Tap Sample ID Prefix field to start editing.
  - Use the on-screen keypad to enter lot file name.
  - Tap < to save new name.
  - Scroll to select Prefix value.
- 6. Tap 🔾 (Home key) to reenter Measurement screen.
- 8. See option to track and log notes.
- 9. From Measurement screen, tap symbol to start logging, and tap symbol to stop logging.

# 12.3. AUTOHOLD LOGGING

- 1. From Measurement screen, tap 🗢 (Measurement Menu).
- Tap Reading tab then select stability criteria (Accurate, Medium, or Fast). Note: Autohold logging uses this criteria for logging. Setting this will affect when data is recorded.
- 3. Tap to select Direct/Autohold reading mode.
- 4. Tap Logging tab and select Autohold logging type.

12:42:49 03/04/2022		Ŕ	<b>•</b> •	0%	HI6321
	Measurement Settings				
Calibration	Logging Type	Automatic	Manual	Autohold	
Reading Temperature	Sampling Period	180 min 1 sec 2 sec			
View	File Name	Create	UnnamedLog		
Alarms	Log Note	Lab 112			
Logging	Log Info 1	test			
Profiles	Log Info 2				

- 5. Create a file name or have a file previously set up.
- 6. Option to define sample ID (name and prefix)
  - Tap Sample ID Prefix field to start editing.
  - Use the on-screen keypad to enter lot file name.
  - Tap < to save new name.
  - Next, scroll to select Prefix value.
- 7. Tap 🔾 (Home key) to reenter Measurement screen.
- From Measurement screen, tap 
   Oirect logging symbol) to start logging. Active Autohold logging in progress is indicated by displayed 
   symbol. Progress bar, top left corner, indicates progress status.
- 9. Once configured stability criteria is checked, and conditions are met, data is logged. The instrument prompts measurement screen, displayed in configured view.

### CSV file format naming convention

To help identify measurement-specific data, files within a data package have the measurement unit automatically included in the file name (e.g. when reading and logging in Conductivity Parameter, automatically logged files retain the "EC" ID in the file name).

Manual or Autohold logged files retain the distinction between different reading modes in the file name. Users have the option to input file name:

- EC > 20220329\_085101-ec\_auto
- Resistivity > 20220309\_095704-res\_auto
- Salinity > 20220331\_101507-**sal**\_auto
- TDS > 20220331\_101458-tds\_auto

# **13. MAINTENANCE**

### **Probe Maintenance**

Rinse the probe thoroughly as water residue may not be visible.

### Cleaning

Dirty or improperly cleaned probes can result in erratic and inaccurate readings.

- Clean off the external sheath with a soft cloth and surfactant solution.
- Rinse the probe under a stream of running tap water to remove salt or minerals. Jet the tap water stream through the opening to dislodge any debris.
- Only if strictly necessary, carefully remove the outer plastic sheath to disassemble the probe. Clean off with a warm water (surfactant) mixture and follow with a through rinsing with purified water. Allow pieces to dry and reassemble.

Calibrate the probe with the appropriate standard solution for the intended application.

### Storage

- Store the probe dry, after cleaning in distilled water.
- Clean the probe and calibrate after long-term storage.

### General Cleaning of HI6321

The following steps outline the process to ensure users keep the benchtop clean and disinfected while limiting the risk of damage from unsuitable cleaners.

- Disinfect the screen using commercially available non-ammonia glass or disinfectant cleaner.
- Apply a small amount of cleaner directly to a microfiber or lint-free disposable cloth. Make sure the cloth is damp and not wet.
- Wipe the glass touch screen clean with the cloth. Do not apply cleaner directly to the interface.

# **14. ERROR MESSAGES**

The instrument shows clear warning messages (bottom of the screen) when erroneous conditions appear or when measured values are outside the expected range.

The information below provides an explanation of the errors and warnings, and recommended action to be taken.

Displayed Message	Explanation & Recommended Action
"Temperature under/over range"	Temperature outside specified range. Replace probe if necessary.
"Under/over compensation range"	During conductivity calibration, the temperature is under/over the conductivity calibration solution compensation limit.
"EC under/over range"	Outside range in the corresponding scale. Ensure the vent holes are completely submerged and the sample is within specified range.
"Broken electrode"	The meter fails to calibrate or gives faulty readings. Replace the probe.
"EC out of calibration range"	Displayed when the measured value is outside calibration range.
"Factory calibration expired"	Contact the Hanna technical support for the periodic factory calibration.

# **15. SOFTWARE UPDATE**

To introduce new features and/or performance improvements Hanna Instruments<sup>®</sup> releases updated firmware versions. To check for new releases, scan the QR code or go to: <u>https://software.hannainst.com</u>.

### Requirements

- USB-A drive (FAT32 format)
- Device must be turned off

#### Steps

- 1. Scroll down the software downloads page to find the Instrument Firmware list.
- 2. Connect the USB-A drive to PC.
- 3. Find the Firmware Version needed for download, then click DOWNLOAD NOW.
- 4. Wait for \*.hiup file download to complete.
- 5. Plug the flash drive into the USB-A port and turn the meter on.



6. Wait for update to complete (~1 hour). *Note:* Do not turn off the meter or disconnect from power.





# **16. ACCESSORIES**

#### **CONDUCTIVITY SOLUTIONS**

HI7030M or HI7030L	12880 $\mu$ S/cm standard solution, 250 or 500 mL
HI7031M or HI7031L	1413 µS/cm standard solution, 230 or 500 mL
HI7033M or HI7033L	84 $\mu$ S/cm standard solution, 230 or 500 mL
HI7034M or HI7034L	80000 $\mu$ S/cm standard solution, 250 or 500 mL
HI7035M or HI7035L	111800 $\mu$ S/cm standard solution, 230 or 500 mL
HI7037M or HI7037L	100 % NaCl seawater salinity standard solution, 250 or 500 mL
HI7039M or HI7039L	5000 $\mu$ S/cm standard solution, 250 or 500 mL
OTHER ACCESSORIES	
HI740036P	100 mL beaker (10 pcs.)
HI764060	Electrode holder
HI900946	115 Vac to24 Vdc power adapter, US plug
HI900947	230 Vac to 24 Vdc power adapter, European plug
HI920016	USB type A to C cable

# CERTIFICATION

All Hanna® instruments conform to the CE European Directives and UK standards.



**Disposal of Electrical & Electronic Equipment.** The product should not be treated as household waste. Instead hand it over to the appropriate collection point for the recycling of electrical and electronic equipment which will conserve natural resources.

Ensuring proper product disposal prevents potential negative consequences for the environment and human health. For more information, contact your city, your local household waste disposal service, or the place of purchase.

# **RECOMMENDATIONS FOR USERS**

Before using this product, make sure it is entirely suitable for your specific application and for the environment in which it is used. Any variation introduced by the user to the supplied equipment may degrade the meter's performance. For yours and the meter's safety do not use or store the meter in hazardous environments.

# WARRANTY

HI6321 is warranted for two years against defects in workmanship and materials when used for its intended purpose and maintained according to instructions. Electrodes and probes are warranted for a period of six months. This warranty is limited to repair or replacement free of charge. Damage due to accidents, misuse, tampering or lack of prescribed maintenance is not covered.

If service is required, contact your local Hanna Instruments office. If under warranty, report the model number, date of purchase, serial number (see engraved on the bottom of the meter) and the nature of the problem. If the repair is not covered by the warranty, you will be notified of the charges incurred. If the meter is to be returned to Hanna Instruments, first obtain a Returned Goods Authorization number from the Technical Service department and then send it with shipping costs prepaid. When shipping any meter, make sure it is properly packed for complete protection.

# **REGULATORY NOTICES FOR THE WI-FI MODULE**

United States (FCC) FCC ID: 2ADHKATWINC1500.

This device complies with Part 15 of the FCC Rules. Operation is subject to the following two conditions: (1) this device may not cause harmful interference, and (2) this device must accept any interference received, including interference that may cause undesired operation.

Canada (ISED) IC: 20266-WINC1500PB HVIN: ATWINC1500-MR210PB PMN: ATWINC1500-MR210PB

This device complies with Industry Canada's license exempt RSS standard(s). Operation is subject to the following two conditions: (1) This device may not cause interference, and (2) This device must accept any interference, including interference that may cause undesired operation of the device. Le présent appareil est conforme aux CNR d'Industrie Canada applicables aux appareils radio exempts de licence. L'exploitation est autorisée aux deux conditions suivantes: (1) l'appareil ne doit pas produire de brouillage, et (2) l'utilisateur de l'appareil doit accepter tout brouillage radioélectrique subi, même si le brouillage est susceptible d'en compromettre le fonctionnement.

Under Industry Canada regulations, this radio transmitter may only operate using an antenna of a type and maximum (or lesser) gain approved for the transmitter by Industry Canada. To reduce potential radio interference to other users, the antenna type and its gain should be so chosen that the equivalent isotropically radiated power (e.i.r.p.) is not more than that necessary for successful communication. Conformément à la réglementation d'Industrie Canada, le présent émetteur radio peut fonctionner avec une antenne d'un type et d'un gain maximal (ou inférieur) approuvé pour l'émetteur par Industrie Canada. Dans le but de réduire les risques de brouillage radioélectrique à l'intention des autres utilisateurs, il faut choisir le type d'antenne et son gain de sorte que la puissance isotrope rayonnée équivalente (p.i.r.e.) ne dépasse pas l'intensité nécessaire à l'établisse-ment d'une communication satisfaisante.

Japan (MIC) 005-101762

South Korea (KCC) R-CRM-mcp-WINC1510MR210P

Taiwan (NCC) CCAN18LP0321T2

注意!依據 低功率電波輻射性電機管理辦法 第十二條 經型式認證合格之低 功率射頻電機, 非經許 可, 公司、商號或使用者均不得擅自變更頻率、加大功 率或 變更原設計 之特性及功能。第十四條 低功率射頻電機之使用不得影響 飛航安全及 干擾合法通信; 經發現有干擾現象時, 應立即停用, 並改善至無干 擾時 方得繼續使用。前項合法通信, 指依電信規定作業之無線電信。低功率射 頻電機須忍受合法通信或工業、科學及醫療用 電波輻射性 電機設備之干擾。

China (SRRC) CMIIT ID: 2018DJ1305

ANATEL 08497-18-08759

Note: FCC information is marked on the back of the device.